

試藥滾動混合器 Super Roller K.L.-600A

操作說明與維修手冊 Easy instructions

1. 儀器部位指引 Instrument parts

- | | | |
|---|-----------|---------------------|
| A | 開關 | Power switch |
| B | 旋轉筒 (6支) | Roller tubes (6 ea) |
| C | 鋁製護牆 | Wall |
| D | 放置腳墊 (4個) | The base (4 ea) |
| E | 保險絲 | Fuse (0.5A) |
| F | 電源線 | Power line |
| G | 調整速度 | Speed adjustment |

2. 使用說明 INSTRUCTIONS

- 開關出現紅燈亮時，表示電源正在使用當中，關機時紅燈是不出現。
Red light indicates device is in operation.
- 速度調控20 ~ 150 (±3) rpm/min.。
Speed regulation between 20~150 (±3) rpm/min.
- 當長時間連續旋轉超過24小時，請讓本儀器休息30min，再行運作。
A 30 minutes pause is recommended for more than 24 hours continuous operation.
- 請在0 ~40℃ 的溫度範圍內進行運轉。
The equipment should be operated at 0~40℃.
- 500mL以下之罐(瓶)直接放在旋轉筒上。需注意罐(瓶)口是否鎖緊，以防液體滲出。
Place tubes or bottles directly onto rolling facilities. Tubes or bottles (under 500 ml) should be tightly closed before exercise to avoid leaking.
- 旋轉筒使用時，切勿用手阻擋旋轉，維持旋轉方向。
When device in operation, do not obstruct cycling.

3. 故障排除 COMMON FAULTS and TROUBLESHOOTING

狀況 Symptoms	可能原因 Check points	排除方式 Shooting
旋轉筒不轉動 (開關即亮紅燈) The switch or red light is on but no rolling.	旋轉筒齒輪受異物卡住或放置超過 3 年未啟用。 可能儀器內部潤滑油能漏出。 Cogwheel being stopped by small objects, for device has not been used for quite a while.	雙手輔助旋轉筒旋轉，如果還是不轉動，請以電話聯絡服務廠商。 Use your hand to push gently. If not working, call our service dept.
旋轉筒不轉動 (啟動開關時，紅燈仍未亮) No power light and not working after device is turned on.	保險絲鬆掉 Socket is loose or faulty fuse. .	更換本機保險絲 check the voltage of socket or replace fuse.

Manufacturer :

 廣林貿易有限公司
WANG LIN SCIENTIFIC TRADING CO., LTD.
高雄海山路111號樓一第34號14樓 (國泰大戲院)
705 No. 34th Floor, 1st 34th Floor, 14th Floor, Gateway
Tel: +886 7 3330485 Fax: +886 7 3330500
http://www.wanglin.com.tw
E-Mail: wanglin@ms26.hinet.net

